



# iCut-40P

## Metallographic Precision Cutting Machine



Video



### Contact us

Mikrosize Precision Instrument Co.,Ltd

A-4035 RuiFeng Business Expo, Wuhu City, China , 241000.

Web:[www.mikrosize.com](http://www.mikrosize.com)

Email: [mikrosize@mikrosize.com](mailto:mikrosize@mikrosize.com)



Web:[www.mikrosize.com](http://www.mikrosize.com)

Email:[mikrosize@mikrosize.com](mailto:mikrosize@mikrosize.com)

# Product Features and Application

## Product Features

- High positioning accuracy;
- Wide speed range;
- Strong cutting capacity;
- Built-in circulating cooling system;
- Preset feed speed;
- Touch screen control display, easy to operate and control;
- Automatic cutting can ensure uniform cutting consistency on the surface of the sample;
- Large and bright cutting guard with safety switch.

## Product Application

### ● Industrial manufacturing field

It can be used for the processing and manufacturing of mechanical parts, and can achieve precise cutting of various flat materials. By precisely controlling the cutting angle and depth, the precision flat plate cutting machine can ensure that the material is cut accurately and completely, meeting the high-precision requirements of industrial manufacturing.

### ● Architectural model industry

Precision flat plate cutting machines can ensure the production needs of architectural models, such as the realization of precise proportions and dimensions.

### ● Medical industry

Precision flat plate cutting machines can produce surgical instruments and equipment that require delicate operations through precise cutting.



**Machine on the Front**



**The Left Side of Machine**

# Instrument Appearance



**1. Transparent protective cover: used to prevent cooling water and cutting waste from splashing, and the transparent material makes it easy to observe the state of the cut sample**

**2. Cooling water valve: used to control the size and switch of the cooling water flow**

**3. Cooling water nozzle**

**4. Clamp**

**5. Closed-door detection**



**1. Start button**


**2. Stop button**

**3. Display screen**

**4. On/off button**

**5. Emergency stop button**

## Instrument Appearance

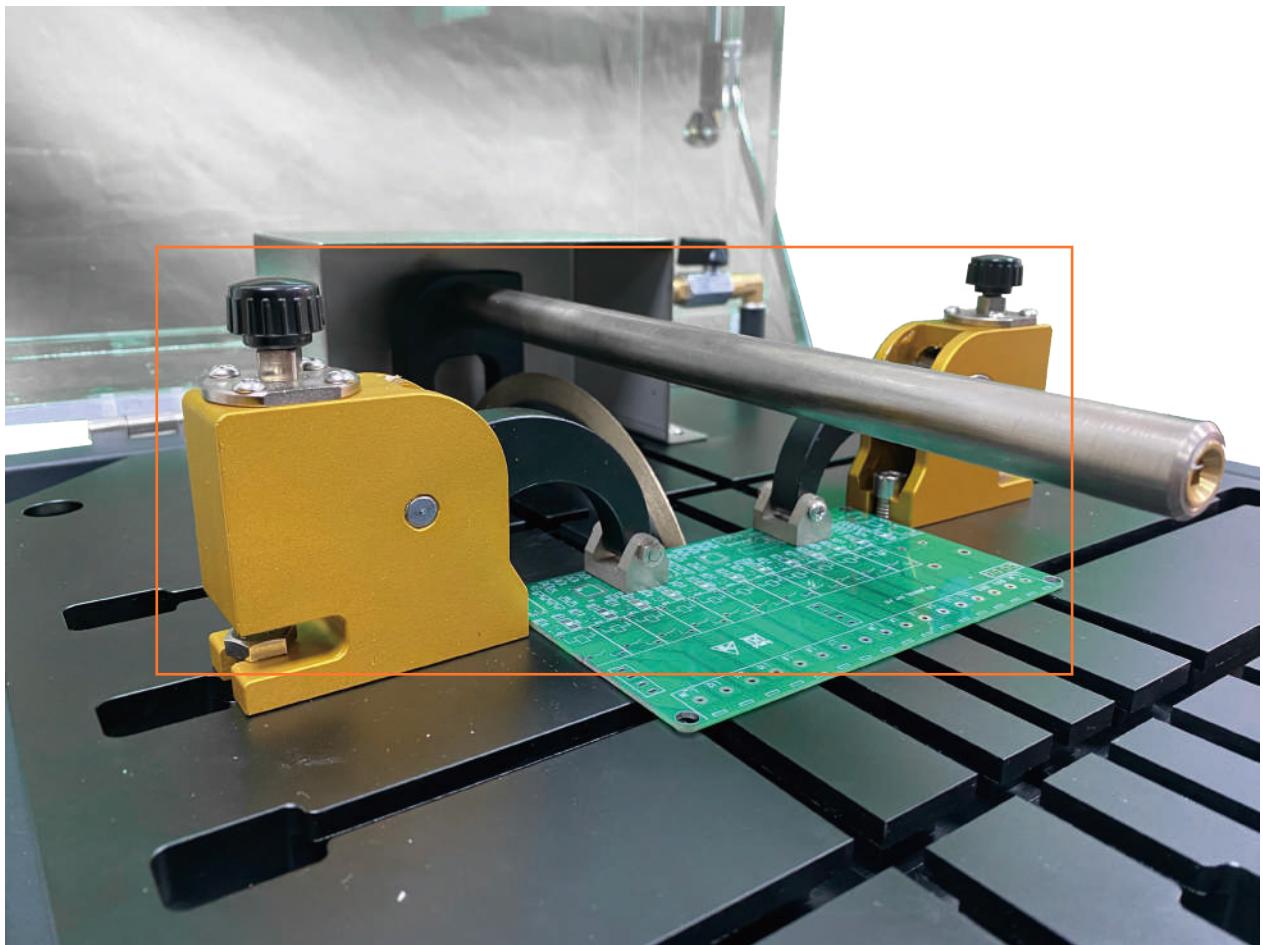

### Clamp

Double-sided quick clamps are used to clamp the specimen quickly and firmly.

# Instrument Appearance



## Cutting Disc


The diamond sintered cutting disc has the advantages of high hardness and high wear resistance.



## Closed-door Detection

Be sure to close the protective cover before starting the equipment. When the protective cover is opened, the system alarm interface will prompt "Door Open". When the equipment is cutting, opening the protective cover will force the machine to stop.

## Instrument Appearance

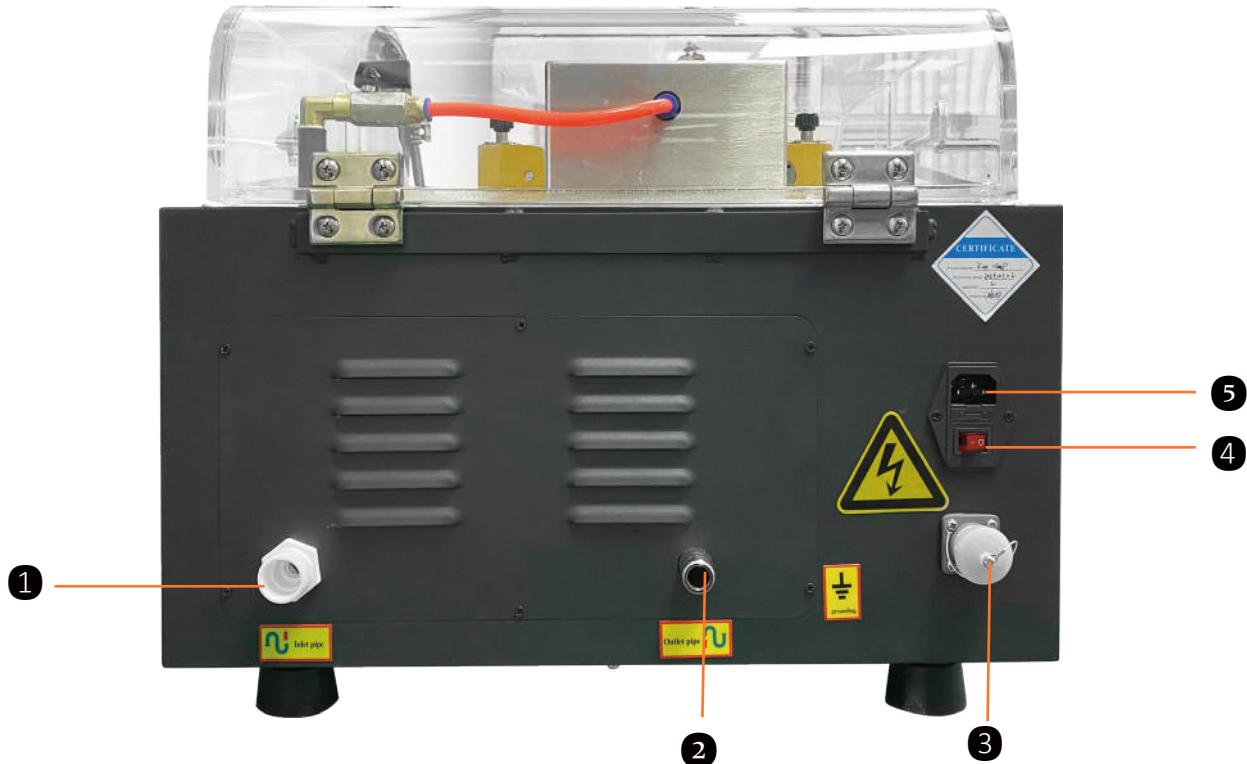

**1. Start Button**

**2. Stop Button**

**3. Display Screen**

**4. On/Off Button**

**5. Emergency Stop Button**

# Instrument Interface & Function

2025-01-13 16:28:55

## CATALOG INTERFACE


WORKING INTERFACE

SETTING INTERFACE

ALARM INTERFACE

ENGLISH

After the device is turned on, it enters the "CATLOG INTERFACE" interface. This device consists of four parts: "WORKING INTERFACE", "SETTING INTERFACE", "ALARM INTERFACE" and "LANGUAGE SWITCHING".

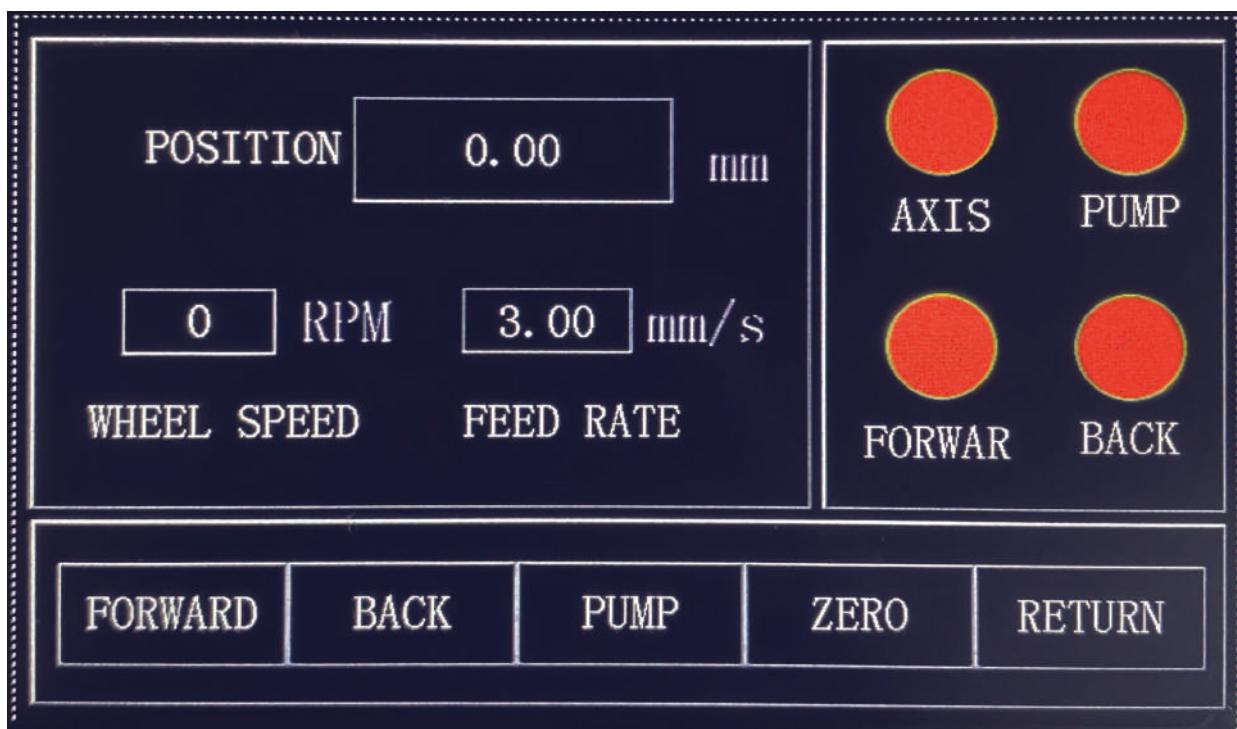

### 1. WORKING INTERFACE

This interface displays the operating status of the equipment, including the status indicators of the spindle, feed axis, water pump, spindle speed and cutting speed.


The FORWARD and BACK buttons drive the feed axis.

The PUMP button turns on the cleaning and cooling water pump.

The ZERO button resets the current position.

The RETURN button returns to the catalog page.

# Instrument Interface & Function

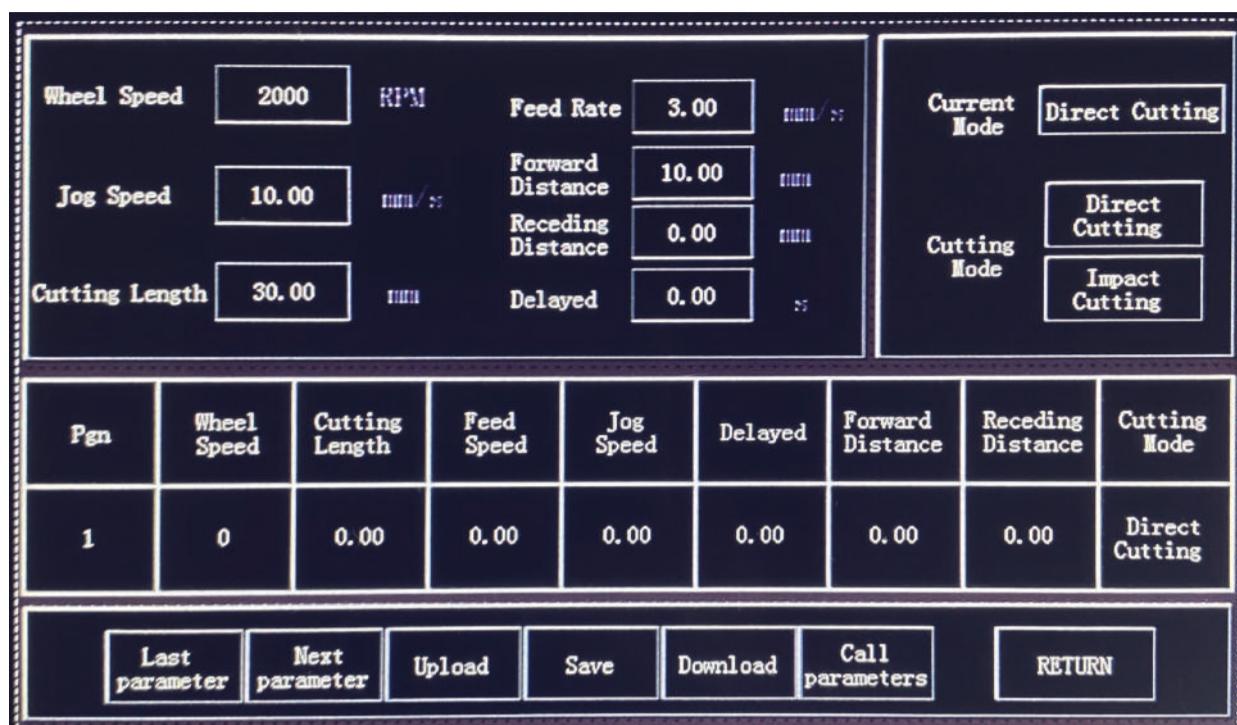

## 2.SETTING INTERFACE

This interface is the parameter setting interface of the device. You can set Wheel Speed, Jog Speed, Cutting Length, Feed Rate, Forward Distance, Receding Distance, and Delayed within the allowed range. Touch the screen to display numeric keys, and enter the required values.

Two cutting modes Direct Cutting and Impact Cutting are available.

The set cutting parameters can be saved and directly called for later use. Up to 10 sets of cutting data can be saved.

Alarm Time	No.	Alarm matt
2025-01-13 16:23:24	3	The door is open
2025-01-13 16:23:24	2	Rear limit
2025-01-13 16:15:14	3	The door is open
2025-01-13 16:11:57	2	Rear limit

Buttons at the bottom:

- REMOVAL
- RETURN


## 3.ALARM INTERFACE

This interface is a record of the device alarm prompt, including the time and content of the alarm.

## 4. LANGUAGE SWITCHING

This device supports both Chinese and English. Touch the screen button to switch the language.


## Instrument Dimension

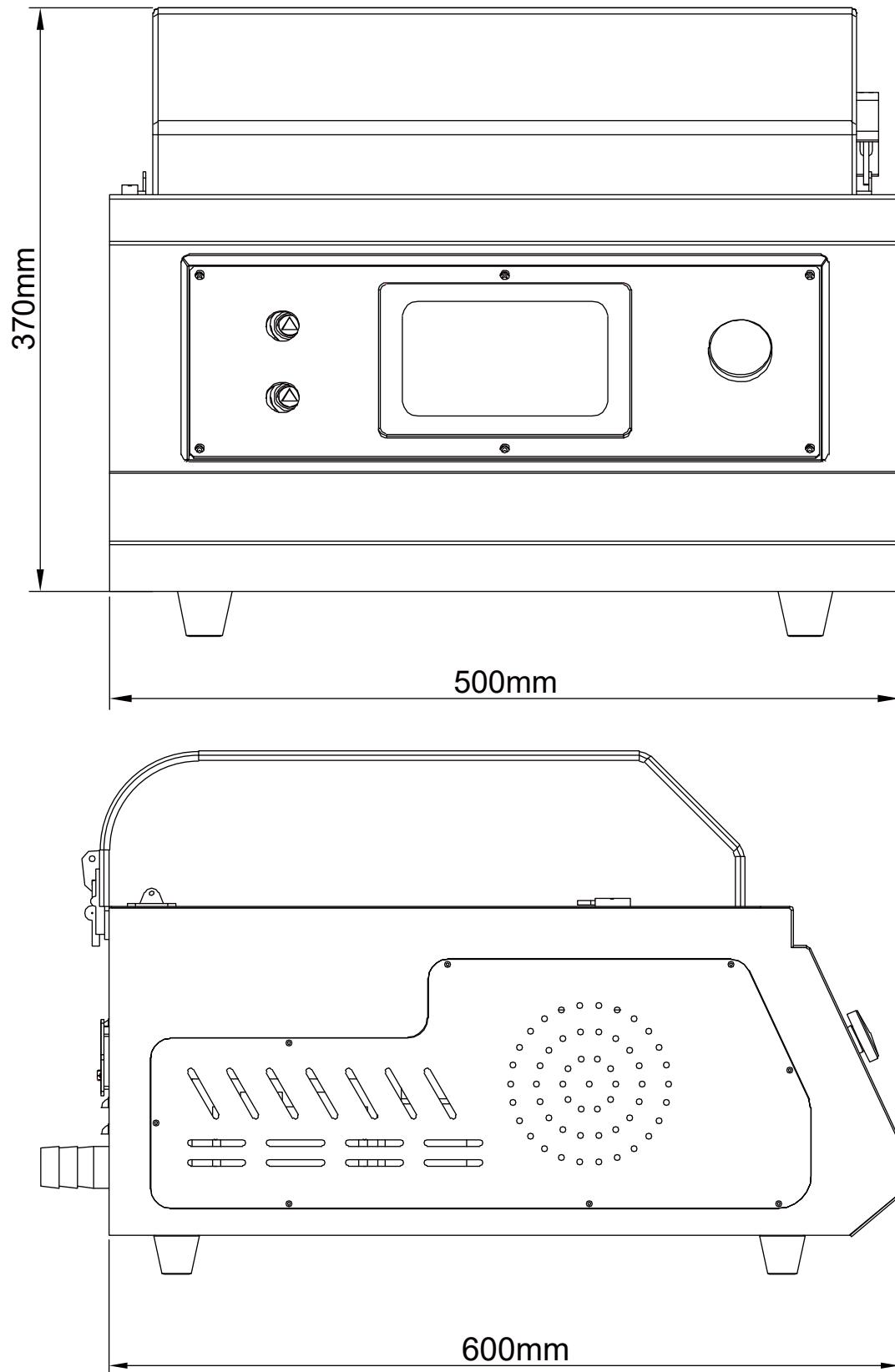

# Technical Specification

<b>Model</b>	iCut-40P
<b>Product Name</b>	Metallographic Precision Cutting Machine
<b>Y-axis Movement</b>	200mm
<b>Cutting Method</b>	Straight Line, Pulse
<b>Cut-off Wheel</b>	φ200*0.9*32mm
<b>Rotation Speed</b>	500-3000r/min, Customized
<b>Automatic Cutting Speed</b>	0.01-3mm/s
<b>Manual Speed</b>	0.01-15mm/s
<b>Impact Cutting Distance</b>	0.1-2mm/s
<b>Maximum Cutting Thickness</b>	40mm
<b>Maximum Clamping Length of Table</b>	425mm
<b>Maximum Clamping Width of Table</b>	200mm
<b>Display</b>	5 Inch Touch All-in-one Control
<b>Data Usage Method</b>	10 Kinds Can Be Selected
<b>Working Table Dimension</b>	425*435mm
<b>Motor Power</b>	600W
<b>Power Supply</b>	Single Phase 220V
<b>Machine Dimension</b>	530*600*410mm
<b>Net Weight</b>	85KG
<b>Gross Weight</b>	101KG

# Standard Delivery

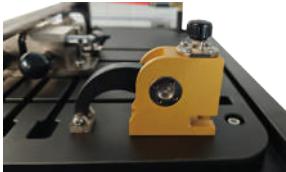Name	Specification	Qty	Picture
Machine Mainframe	iCut-40P	1 set	
Cut-off Wheel	φ200*0.9*32mm	1 pc	
Inlet Pipe		1 pc	
Outlet Pipe		1 pc	
Clamp	Choose 1 clamp in the standard delivery	1 pc	 
Wrench	17-19mm	1 pc	
Hex Wrench	M4	1 pc	
Instruction Manual		1 copy	
Warranty Card		1 copy	
Product Certificate		1 copy	